




GC CAPILLARY COLUMN APPLICATIONS: Industrial Solvents

C1 to C10 Ethyl Esters

COLUMN: BTR-CW, Bonded Carbowax/BTR - Oxygen Resistant 30M. x 0.25mm I.D. x 0.25µm film
Cat. No.: BTR-CW-30-0.25F

Temperature: 80° (4 min. hold) (8°/min.) - 200°C
Injector: 250°C
Detector: 250°C
Split Flow: 40cc/min.
Sample: 0.4µl

- | | | |
|-------------------|-------------------|---------------------|
| 1. C1 ethyl ester | 5. C4 ethyl ester | 9. C8 ethyl ester |
| 2. C2 ethyl ester | 6. C5 ethyl ester | 10. C9 ethyl ester |
| 3. C3 ethyl ester | 7. C6 ethyl ester | 11. C10 ethyl ester |
| 4. solvent | 8. C7 ethyl ester | |

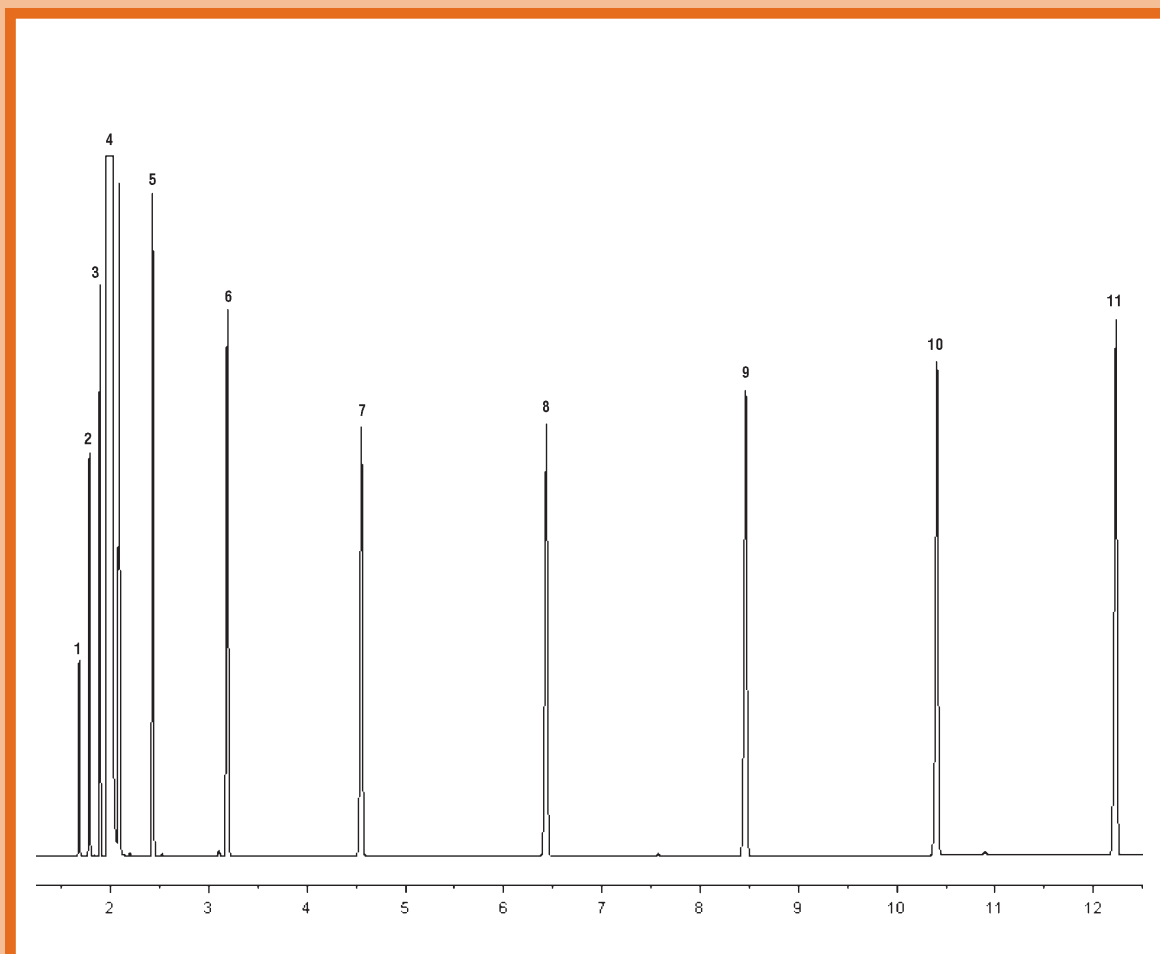

QUADREX



PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com • WEB: www.quadrexcorp.com