



GC CAPILLARY COLUMN APPLICATIONS: **Clinical/ Toxicological**


Mace

Column: 007-17, (50% Phenyl) Methylpolysiloxane
25M. x 0.25mm I.D. x 0.25µm film
Cat. No.: 007-17-25-0.25F

Temperature: 100° (1 min. hold) (10°/min.) -280°C
Injector: 220°C
Detector: MSD
Carrier Gas: 28 cm/sec., Helium

1. 2-Chloroacetophenone
2. contaminant

3. contaminant
4. o-Chlorobenzalmononitrile

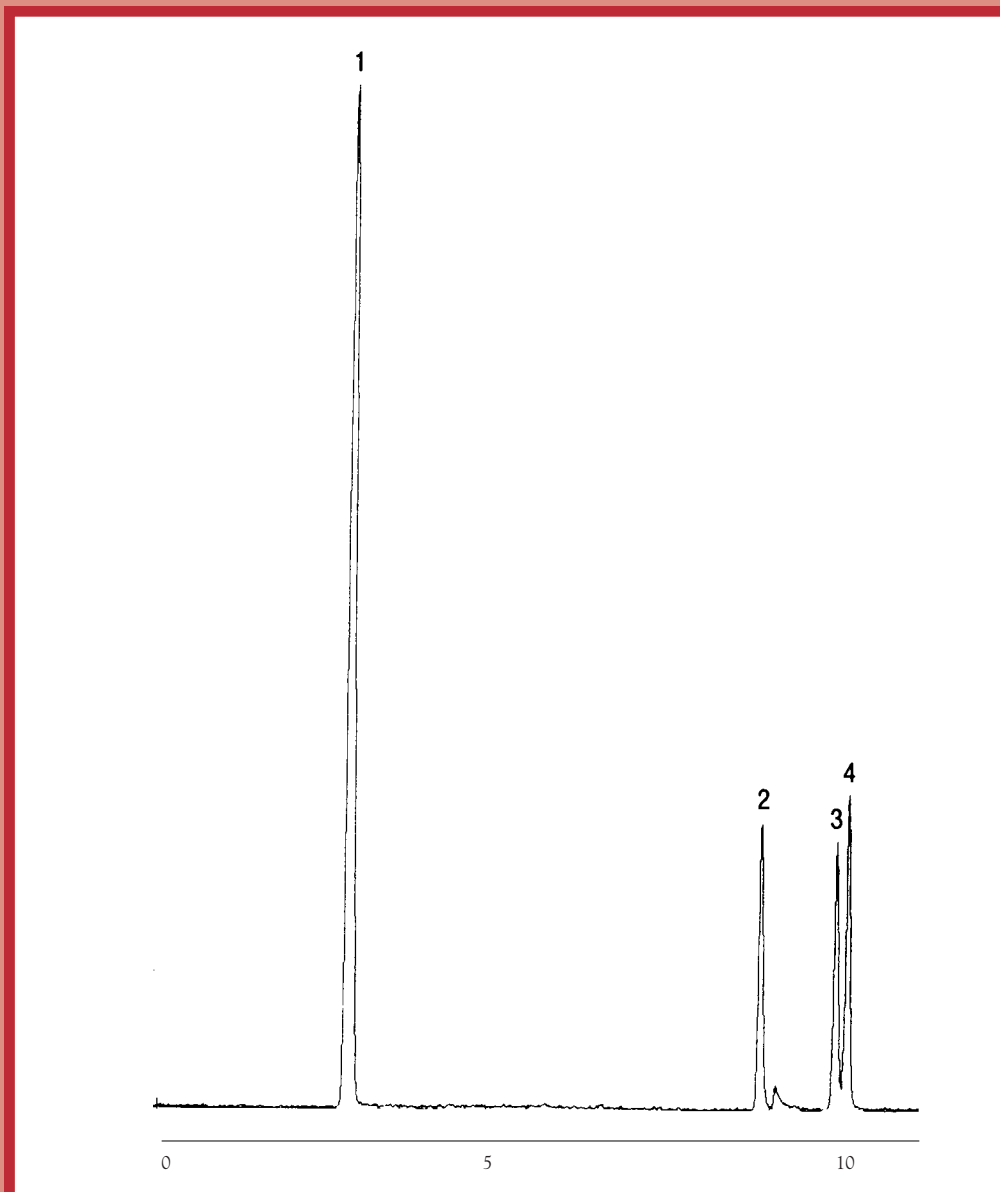

QUADREX



PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com • WEB: www.quadrexcorp.com