



# GC CAPILLARY COLUMN


## APPLICATIONS: Ultra Alloy™ Stainless Steel

### *Residue Solvents in Paint*

**COLUMN:** UAC-CW, Carbowax  
30M. x 0.50mm I.D. x 1.0µm film  
Cat. No.: UAC-CW-30V-1.0F

**Temperature:** 50° (3 min. hold) (5°/ min.) - 100°C  
**Injector:** 200°C  
**Detector:** 200°C, FID  
**Carrier Gas:** 3.4 ml/min., Helium

1. THF                      3. MEK                      5. toluene  
2. ethylacetate        4. iso-C3H7OH

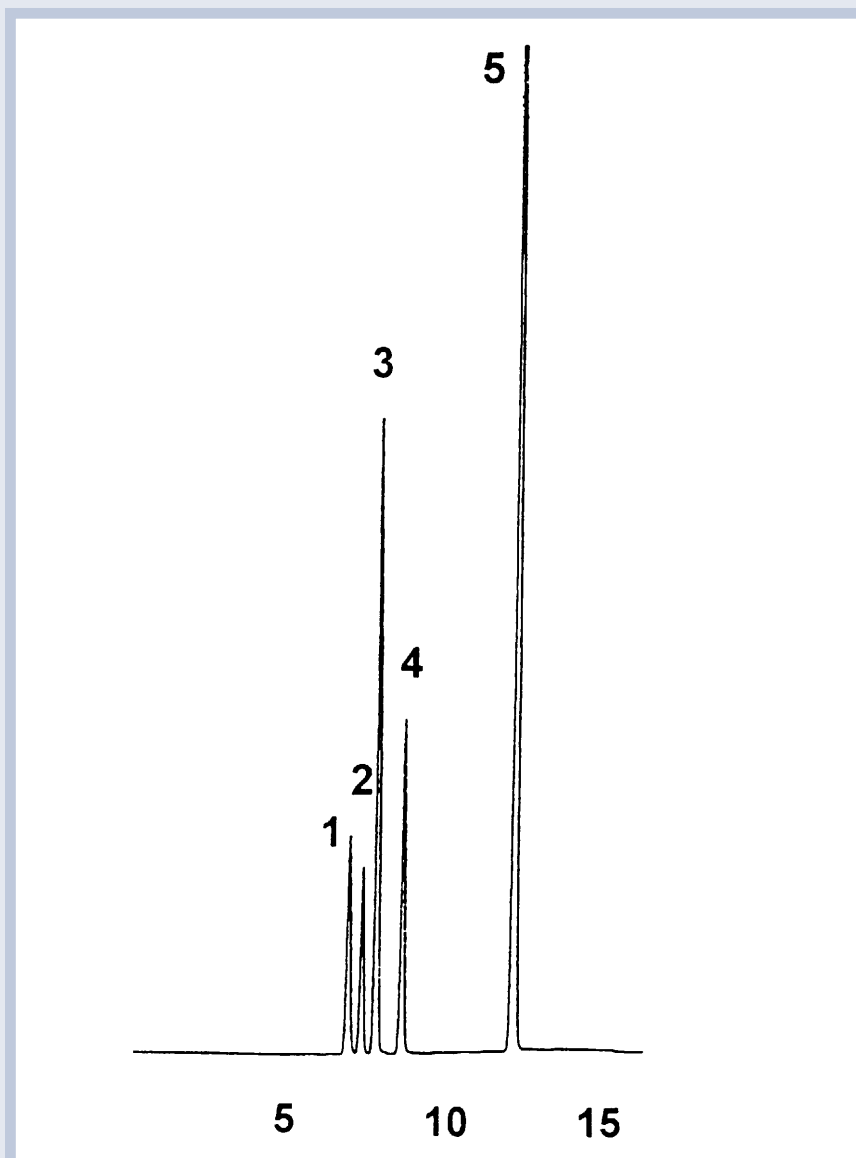

**QUADREX**



PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391  
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com • WEB: www.quadrexcorp.com