



GC CAPILLARY COLUMN

APPLICATIONS: Ultra Alloy™ Stainless Steel

Solvent Mixture

COLUMN:

UAC-1, Dimethylpolysiloxane
60M. x 0.25mm I.D. x 1.0µm film
Cat. No.: UAC-1-60-1.0F

Temperature:

50° (15 min. hold) (10°/min.) - 250°C

Injector:

200°C


Detector:

300°C

1. methanol
2. ethanol
3. acetone
4. isopropyl alcohol
5. dichloromethane

6. ethylacetate
7. chloroform
8. tetrahydrofuran
9. n-butyl alcohol
10. pyridine

11. dimethylformamide
12. cyclohexanone
13. phenethylamine

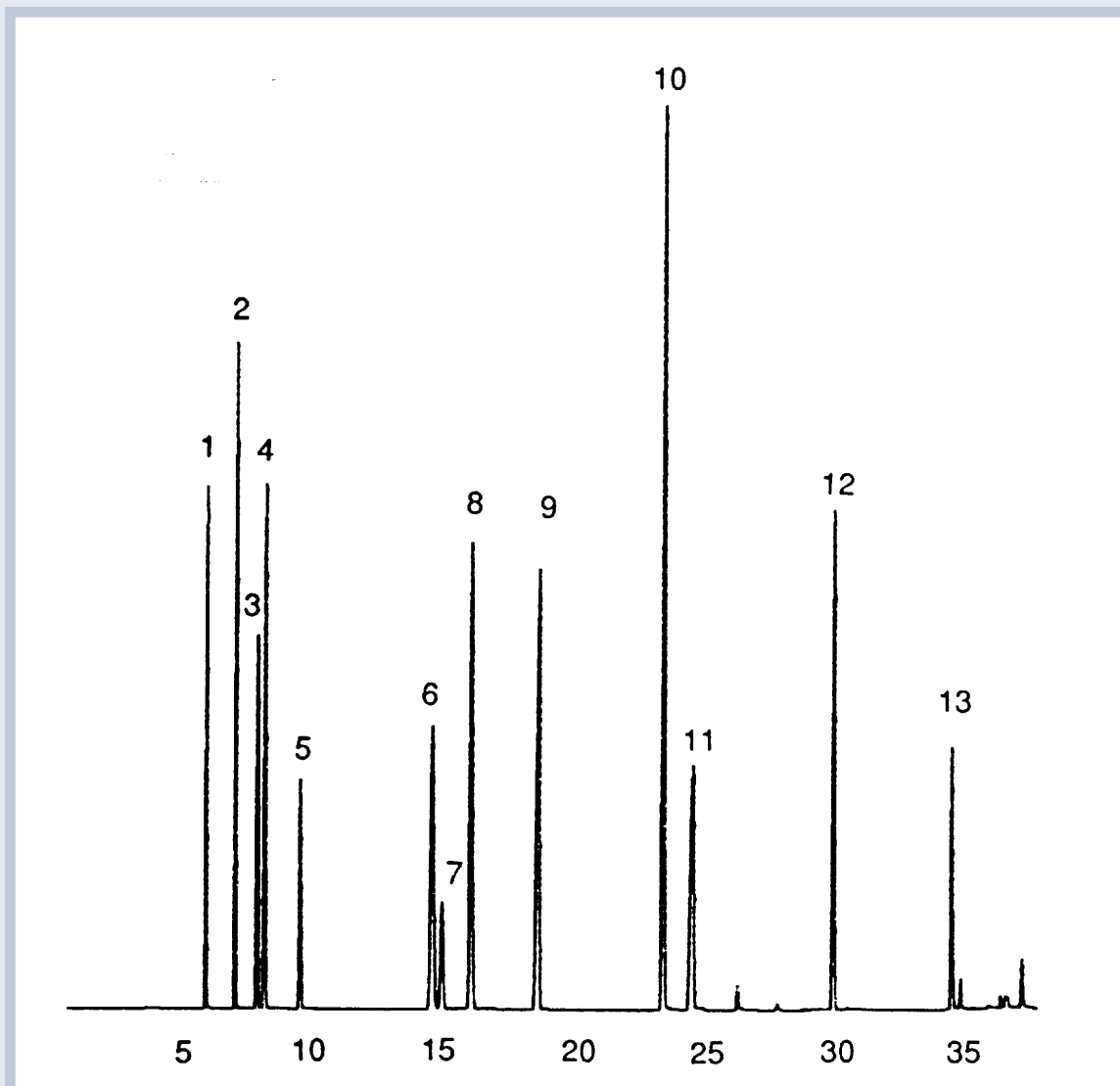

QUADREX



PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com • WEB: www.quadrexcorp.com