




GC CAPILLARY COLUMN APPLICATIONS: Fames

*Volatile
Free
Fatty
Acids*

COLUMN: 007-FFAP, Bonded FFAP
15M. x 0.32mm I.D. x 0.25µm film
Cat. No.: 007-FFAP-15W-0.25F

Temperature: 110° (10°/min.) - 250°C
Injector: 220°C
Detector: 250°C, FID
Carrier Gas: 26 cm/sec., Helium

- | | | | |
|---------------|---------------|--------------|---------------|
| 1. acetic | 5. valeric | 9. caprylic | 13. palmitic |
| 2. propionic | 6. isocaproic | 10. capric | 14. stearic |
| 3. butyric | 7. caproic | 11. lauric | 15. arachidic |
| 4. isovaleric | 8. enanthic | 12. myristic | |

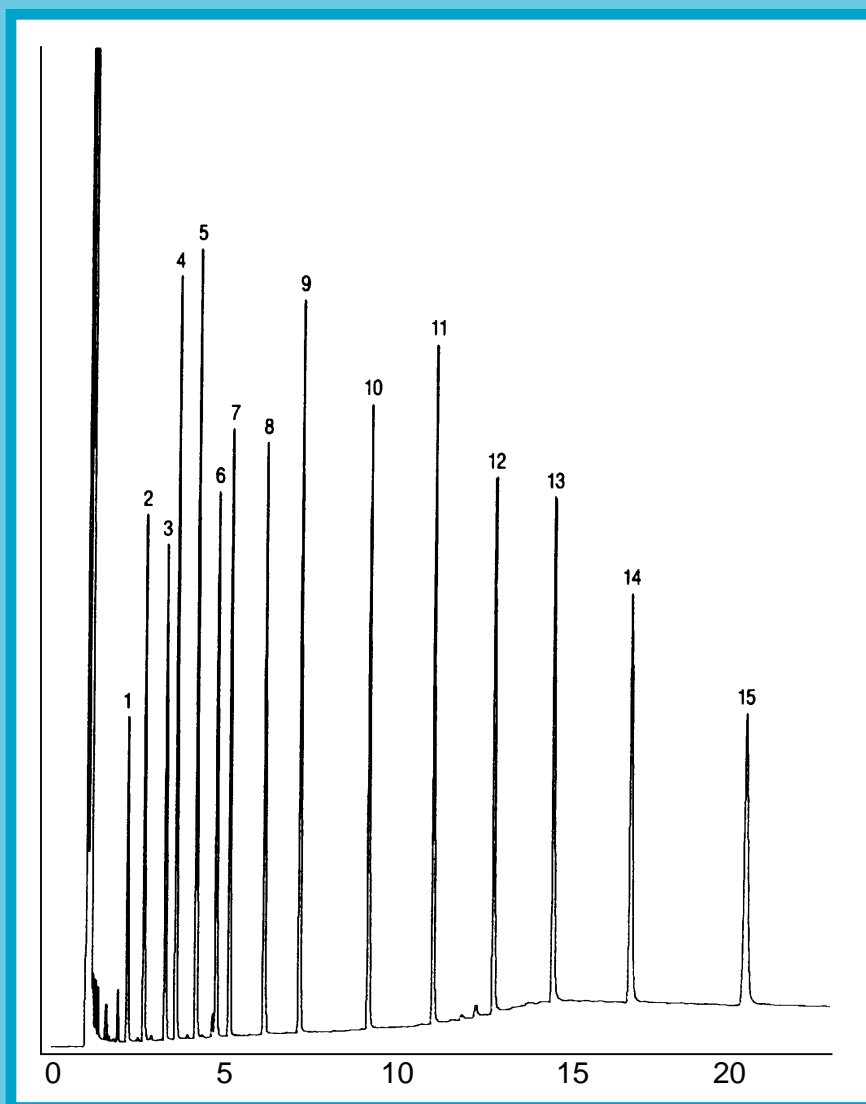

QUADREX



PO Box 3881, Woodbridge, CT 06525 • (203) 393-3112 • FAX (203) 393-0391
TOLLFREE: 800-275-7033 • E-MAIL: tech@quadrexcorp.com • WEB: www.quadrexcorp.com